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Polymer-based Aqueous SEC (GFC) Columns for Bionanoproduct Analysis and Preparative Use: UB series

UB NEW	Suitable pore design for separating bionanoproducts such as viruses, virus-like particles (VLPs), and exosomes
---	--

(Standard columns) [Housing material: SUS]

Product code	Product name	Plate number (TPN/column)	Particle size (μm)	Target Molecular Size (nm)	Column size (mm) I.D. x Length
F6429301	UB-50 NEW	$\geq 3,500$	26	< 50	8.0 x 300
F6429302	UB-100 NEW	$\geq 3,500$	27	> 50	8.0 x 300

Product name	Maximum pressure (MPa/Column)	Flow rate (mL/min)		Temperature range ($^{\circ}\text{C}$)	pH range	Shipping solvent
		Recommended	Maximum			
UB-50, UB-100	0.5	0.5 - 1.0	2.0	4 - 60	2 - 14	H ₂ O

Eluent conditions

- Buffers and aqueous solutions of different salts can be used separately or together. Usable buffers include phosphate, acetate, citrate, and Tris. Usable aqueous salt solutions include sodium chloride, sodium sulfate, potassium sulfate, and ammonium sulfate. Their recommended concentration ranges are 0.05 to 0.3 M.
- Up to 50 % (v/v) methanol, ethanol, isopropanol and acetonitrile are usable.
- Aqueous solutions of urea, guanidine hydrochloride, and dithiothreitol (DTT) which are the protein denaturants can be used up to 8 M, 6 M, and 5 mM, respectively. However, their required concentration are usually high, which may damage the column during the solvent replacement. Thus, it is recommended to dedicate the column for these specific uses.
- SDS surfactant can also be added up to 2% (w/v). However, since surfactants tend to remain on the column, solvent replacement after their use takes a longer time than replacement from general solvents. The replacement time can be shorten by using 30 to 50 % (v/v) methanol.

[Operation Manual / Certificate of Analysis](#)

Operation Manuals and Certificate of Analysis / Inspection Certificate for the following products can be downloaded here.

UB-50

UB-100

[Product Name Index](#)

UB-50

UB-100

[Related Products](#)

Silica-based Aqueous SEC (GFC) Columns: PROTEIN KW-800

High Performance Semi-micro Aqueous SEC (GFC) Columns: KW400

Silica-based Aqueous SEC (GFC) Column for Analysis of Antibody Protein: PROTEIN LW

Polymer-based Aqueous SEC (GFC) Columns: OHpak SB-800 HQ and SB-2000

MALS Compatible Columns for Aqueous SEC (GFC): OHpak LB-800

[Product Information \(Related Information\)](#)

Selection of Aqueous SEC (GFC) Columns

Eluent Conditions for PROTEIN KW Series and LW Series

Eluent Conditions for OHpak SB-800 HQ and LB-800

Features of LB-800 Series (Comparison of Baseline Noise with Using MALS Detector)

Separation Features of UB Series

Durability Against Alkaline Solvents (UB Series)

[General Information on Packed Columns](#)

introduction of column cleaning procedure and column test conditions etc.

Packing Materials for Shodex Packed Columns

General Handling Precautions for Packed Columns

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